

FormaSet™ is a proprietary granular used for vapor control and waste management of spilled aldehydes including formaldehyde, gluteraldehyde and formalin.

Safe and effective, FormaSet neutralizes and absorbs hazardous aldehyde spills by first suppressing vaporization followed by the aldehyde changing into a polymeric compound for safe and easy clean up.

The FormaSet spill kit contains all of the basic personal safety wear and clean-up equipment for quick response and ease of use.



BENEFITS

- Neutralizes and absorbs hazardous aldehyde spills
- Designed for quick response and disposal
- Complies with OSHA standards
- Safety risks and environmental hazards minimized

FEATURES

- FormaSet™ Formaldehyde/ Gluteraldehyde spill powder
- Easy-grip, easy-pour bottle
- A complete kit that includes safety wear and clean-up equipment

OTHER

- FormaSet is also available in pillow, sock, or bulk form

APPLICATIONS

Because formaldehyde vapor is a suspected carcinogen, FormaSet was specifically developed to neutralize and absorb Formaldehyde, Gluteraldehyde and Formalin spills as quickly and as safely as possible.

Following an incident, simply apply the FormaSet granular until the spill is rendered into a dry state. For added safety, a vapor barrier sheet (supplied with the spill kit) should be placed over the spill and remain in place for a minimum of one hour. If a formaldehyde odor continues to be present, additional FormaSet™ should be applied. (SorbaSet S™ can also be applied if increased thickening is desired).

Once in a polymeric form, the resultant product is no longer hazardous and can be removed and easily disposed of using the clean-up equipment supplied with the spill kit.

REGULATORY CONSIDERATIONS

Personnel safety risks and environmental hazards are minimized.

Handles "Foreseeable Emergency during uncontrolled release of hazardous material."
OSHA 1910.1200.

Satisfies spill neutralizer requirements in OSHA 1926.441 & 1910.178. "Facilities must have means to neutralize spilled hazardous materials."

EPA: "New Technology and Spill Containment Programs" Federal Register Vol 55, No. 72/
April 13, 1990; "Agents are available that solidify, neutralize, and suppress vaporization.")



Corporate Offices

P: 603.382.8481
F: 603.378.0816

Pennsylvania Offices

P: 570.848.4186
M: 570.371.8464

www.sorb-tech.com

info@sorb-tech.com

SORBTECH Mfg, Inc.